NMR data processing software

Delta

NMR Software

v5.0

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①

②

Backup Name	Start	End	Status
backup	July 23, 2019 17:37:11	July 23, 2019 17:37:11	Idle (Next Start : July 23, 2019 17:38:11)

Configuration

General Schedule Sub Folders Options

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Remote Maintenance Tool Version 1.1.3 © 2009 - 2017 JEOL RESONANCE

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The screenshot displays the 'Backup Management' interface. At the top, there is a blue header bar with the text 'Backup Management'. Below this is a section titled 'Backup Entries' which contains a table with the following data:

Backup Name	Start	End	Status
backup	July 23, 2019 15:41:08	July 23, 2019 15:41:08	Idle (Next Start : July 23, 2019 15:42:08)

Below the table is a 'Configuration' section with four tabs: 'General', 'Schedule', 'Sub Folders', and 'Options'. The 'Schedule' tab is currently selected. It contains the following settings:

- Schedule Type:** Radio buttons for 'Every Week/Date/Time' and 'Interval'. The 'Interval' option is selected.
- Days:** An empty text input field.
- Hours:** An empty text input field.
- Minutes:** A text input field containing the number '2'. A red arrow points from the number '2' to the right, towards the explanatory text.

At the bottom of the configuration section, there are five buttons: 'Add', 'Modify', 'Pause', 'Resume', and 'Delete'.

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The screenshot displays the 'Backup Management' interface. At the top, there is a 'Backup Entries' table with the following data:

Backup Name	Start	End	Status
backup	July 23, 2019 15:43:08	July 23, 2019 15:43:08	Idle (Next Start : July 23, 2019 15:44:08)

Below the table is the 'Configuration' section with tabs for 'General', 'Schedule', 'Sub Folders', and 'Options'. The 'Options' tab is selected. Under the 'Log' section, there is a 'Log File' input field and a list of options: Log Backup Start/End, Log Copy Start, Log Copy Result, and Log Errors. Under the 'Miscellaneous' section, there is a 'Minutes After' input field containing the value '2', with a red arrow pointing to it from the text '表示新创建的数据在2分钟后会自动备份'. Below this is an 'Options' section with Skip Error Files. At the bottom left, the 'Add' button is highlighted with a red box. At the bottom right, there are buttons for 'Modify', 'Pause', 'Resume', and 'Delete'. The footer text reads: 'Remote Maintenance Tool Version 1.1.3 © 2009 - 2017 JEOL RESONANCE Inc. All rights reserved.'

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